Digital Melting Point Apparatus

2935



Year Warranty

Advantages

- TFT Display of Capillary View with High Resolution Viewing Camera
- Extremely Useful for Pharmaceutical Industry
- Highly Accurate & Easy to Operate
- 4 Digit Seven Segment Red LED Display
- Digital Resolution for Temperature 0.1°C
- Silicon oil-bath and in-built magnetic stirrer for uniform heating

Melting Point Apparatus with TFT display & 4 digit LED display is designed for accurate measurement of melting points and for checking the melting range of the solid sample. The apparatus consists of a silicon oil-bath of checking the melting range of the solid samples. The apparatus consist of silicon oil-bath of about 200 ml capacity, covered with a teflon cover with 5 holes, fitted with a special type of oil immersion heater and having three holes for holding the capillaries, one for boiling point tube and one for the thermometer. It is supplied with high resolution camera & TFT display screen to view sample in capillaries.

The apparatus consists of an in-built magnetic stirrer, an electronics controller with indicator for precisely adjusting the rate of heating of silicon oil-bath an ON/OFF switch, a pilot lamp, a magnifier, glare free background illuminator and a High/Low range temperature selector switch and digital temperature indicator with PT-100 RTD sensor.

The 'Start' switch is used to start the measurement of melting point. User can stop the process of measurement using 'Stop' switch after the instrument reaches the melting point of the sample. The display will always show the last temperature where 'Stop' switch was pressed in last.



Digital Melting Point Apparatus

Technical Data	
TEMPERATURE C	0-300 °C on a 4 digit LED display
DISPLAY	TFT screen with high resolution camera for viewing capillaries & Temperatue in 7 segment 4digit LED display
RESOLUTION	0.1 °C in model 2935,935 or 1°C in model 934
CAPILARY GLASS TUBE	75mm long, 0.8 to 1.1mm internal diameter
BOILING POINT TUBE	Glass Tube 6 x 75mm Long
AMOUNT OF SAMPLE	Fine powder properly dried at a temperature considerably below its melting point, 4 to 6 mm high
OIL-BATH	Silicon oil bath beaker - 200 ml capacity with teflon cover fitted with oil immersion heater. It is having three holes for holding the capillaries, one for boiling point tube and one for thermometer
STIRRER SPEED	Controllabe with front knob
POWER	230V ±10% AC, 50 Hz
DIMENSION	370 x 290 x 290mm (LxBxH) (Approx.)
WEIGHT	12 Kg. (Approx.)
ACCESSORIES	TFT Display Screen & High Resolution Camera Box of Capillaries Glass Tubes for Liquid Sample (6 x 75 mm) - 2Nos. Magnetic Capsule Bottle of Silicon Oil - 250 ml Glass Cylindrical Beaker with Teflon cover fitted with immersion heater Instruction Manual Dust Cover

Specifications are subject to change due to continuous

Product Range

Semi Auto Analyser, Double Beam UV/VIS Spectrophotometer, Single Beam UV/VIS Spectrophotometer, Flame Photometer, Heamoglobin Meter, tal Photo Colorimeter, pH Meter, Conductivity Meter, TDS Meter, DO Meter, Salinity Meter, Turbidity Meter, Nephelometer, Colony Counter, Karl Fischer Moisture Titrator, Photo Fluorometer, Dissolution Test Apparatus, Disintegration Test Apparatus, Friability Test Apparatus, Melting Point Apparatus, Telethermometer, Temperature Indicator, Water & Soil Analysis Kit, Portable pH Meter, Portable Conductivity Meter, Portable TDS Meter, Portable DO Meter, Student Microscope, Binocular microscope, Co-Axial Binocular, Medical Microscope

